

# Advanced Test Equipment Corp. www.atecorp.com 800-404-ATEC (2832)

Product data sheet



# FP89-HL Ultra-Low Refrigerated-Heating Circulator

# Upper middle class with ext. Pt100 sensor connection

Ultra-low Refrigerated Circulator with 2-stage cascade-cooling

JULABO Ultra-Low Refrigerated Circulators for heating and cooling are suitable for external temperature tasks. Powerful circulating pump systems and high heating and cooling capacities guarantee short heat-up and cool-down times. ACC 'Active Cooling Control' provides active cooling control across the whole temperature range. An energy saving feature with only minor loss of heat is provided by the proportional cooling capacity control in the 'FP' refrigeration units. The inevitable dust accumulation that occurs in refrigeration systems is solved by means of an easily removed venting grill, which allows for easy cleaning. The units are equipped with handles or castors for an easy transportation. All models have a drain tap on the front allowing the fluid to be drained easily. Additionally the instruments are equipped with an improved insulation to help avoid ice-formation, and have a visual liquid level display.



### Your advantages

- VFD COMFORT DISPLAY
- LCD DIALOG DISPLAY backlit for convenient interactive operation
- Keypad for setpoints, warning/safety values and menu functions
- ICC (Intelligent Cascade Control), self-optimizing temperature control
- TCF Temperature Control Features to optimize the control behaviour
- ATC3 3-Point-Calibration
- Pt100 External sensor connection for measurement and control
- SMART PUMP, electronically adjustable pump stages
- Early warning system for low liquid level (DBGM 203 06 059.8)
- Adjustable high temperature cut-out, visible via display
- RS232/RS485 interface for online communication
- Integrated programmer for 6 x 60 program steps
- Connections for solenoid valve and HSP booster pump
- · Active Cooling Control

#### **Technical Data**

Order No.	9312689
Model series	HighTech
Category	Ultra-Low Refrigerated-Heating Circulators
Working temperature range (°C)	-90 100
Temperature stability (°C)	±0.02

JULABO USA Inc. | 884 Marcon Boulevard | Allentown, PA 18109

Page 1 of 4

### Product data sheet



	THE TEMPERATURE DIRECTOR DISCOUNTY
Setting / display resolution	0.01 °C
Integrated programmer	6x60 steps
Temperature Display	VFD LCD
Heating capacity (kW)	1.3
Cooling capacity (Medium Ethanol)	
20°C	1 kW kW
0°C	0.92 kW kW
-20°C	0.88 kW kW
-40°C	0.75 kW kW
-60°C	0.58 kW kW
-80°C	0.2 kW kW
Pump capacity flow rate (I/min)	22-26
Pump capacity flow pressure (psi)	5.8-10.2
Pump capacity flow suction (psi)	2.9-5.8
Bath opening / bath depth (W x L / D inch)	5.1 x 5.9 / 6.3
Pump connections	M16x1
Barbed fittings diameter (inner dia. / mm)	8 / 12
Filling volume liters	8
Refrigerant	R404AR508B
External Pt100 sensor connection	integrated
Digital interface	RS232 RS485 Optional Profibus
Ambient temperature	540 °C
Dimensions W x L x H (inch)	21.7 x 23.6 x 36.2
Weight (LBS)	297
Classification according to DIN12876-1	Classification III (FL)according to DIN 12876-1
Included with each unit	2 each barbed fittings for tubing 8 and 12 mm inner dia. (pump connections M16x1 female).
Cooling of compressor	Air

# **Available voltage versions**

230 V / 50 Hz 230 V / 60 Hz



#### **Characteristics**



#### A perfect view

Ample, easy to read VFD Comfort display for simultaneous display of 3 values, warning functions, high temperature cut-off, pump stages (resolution 0.01 °C)



# Additional plain text information

Comfortable LCD dialog display for interactive operation with plain text display



#### Pump stage and liquid level

Backlit indicator for selected pump stages and filling volume on Presto® PLUS, Magnum 91 & Forte HT



#### Comfortable and detailed

Comfortable keypad with additional menu functions for pump stages, calibration, control parameters, programmer, warnings, etc.



#### For perfect results

'Intelligent Cascade Control', automatic & self optimizing adjustment of PID control parameters, temperature stability ±0.005 °C internal, <±0.05 °C external



#### **Full control**

'Temperature Control Features' for individual optimization, access to all important control parameters, additional settings for band limit, limits, Co-Speedfactor etc.



## **Highest measuring accuracy**

'Absolute Temperature Calibration' for manual compensation of a temperature difference, 3-point calibration



### **Clever pump system**

Reliable and consistent pump capacity, electronically adjustable pump stages



# Control from the external application

External Pt100 sensor connection for precise measurement and control directly in the external application



# **Connection compliant to standard**

RS232/RS485 dual-interface for serial data transmission according to EIA-485 industry standard (2-wire bus technology), upgradable with Profibus DP



### **Optimal program control**

For the execution of time and temperature dependant profiles, 6 temperature profiles with 60 steps max., with real time clock



# Automatic control of operating time

Electronic countdown-timer function for timer-programmed unit shut-down, standby mode after programmed time expires



### **Energy saving cooling**

Proportional cooling control for automatic adjustment of cooling power or temporary switch-off of compressor as needed to save up to 90 % energy in comparison to unregulated cooling machines



### Consistent cooling capacity

Easily removable venting grid for quick and easy cleaning

#### Product data sheet





# 100 % Cooling capacity

'Active Cooling Control' for cooling available throughout the entire working temperature range, fast cool-down even at higher temperatures



# **Condensation and ice protection**

A heated cover plate prevents condensation or ice build-up in the bath

#### Other characteristics

Optional: Electronic module with analog connections for external programming, temperature recorder, standby, alarm (Order no. 8 900 100).

Technical changes without prior notification. Images may deviate from the original.